THE SMART SOLUTION

Microscope Set 900

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NOTE: CAUTION ADVISED

Children should work with this instrument only under adult supervision. It contains instruments with needle points and sharp cutting edges, such as a scalpel and a slicer.

Read these instructions carefully before use and follow all the directions described in them. We recommend that you retain these instructions. Small children and animals should not be in the area when experiments are being carried out. Store the microscope in a location out of reach of small children.

INSTRUCTIONS FOR PERSONS PROVIDING SUPERVISION

- a) The instructions, safety rules and first aid information need to be obeyed under all circumstances. They should be carefully read and retained for later consultation.
- b) You should only prepare the specimens described in the instructions.
- c) This microscope set is suitable for children aged 8 and over.
- d) Since the capabilities of children can also differ greatly within an age group, persons providing supervision should pay close attention to which specimens are suitable and safe to use in each case. These operating instructions help adults in assessing whether a specimen is suitable for a particular child.
- e) The persons providing supervision should discuss the safety instructions with the children before the preparation of specimens. Special attention must be paid to the safe handling of pointed and sharp objects such as needles, scalpels and cutting instruments.
- f) During the preparation of specimens, the work surface must be free of objects and food that could hinder work. It should be well lit and located near a water connection. A fixed table with a heat-resistant surface should be provided.
- g) Keep a special rubbish bin or a tin can on hand for solid waste.

SAFETY INSTRUCTIONS

General first aid information

- a) In case of burns: Wash the affected area for five minutes using plenty of water.
- b) In case of cuts: Wash the cut with antiseptic solution or use clean water if no such solution is available. Then cover the wound with a bandage. In case of larger cuts, seek medical assistance. A doctor should always be consulted in case of injuries.

SAFETY MEASURES

- a) Read these operating instructions carefully before use of the instrument and follow all directions. Keep these instructions in a safe place.
- b) Small children and animals should be kept away from the area where experiments are being carried out.
- c) Store the microscope set outside of the reach of small children.
- d) Always clean the instruments after use.
- e) Wash your hands after you prepare specimens.
- f) Never use instruments not provided with this set.
- g) You should not drink, smoke or eat near the area where experiments are carried out.
- h) Do not place food back into the original packaging. Throw it away immediately.


PRECAUTIONARY MEASURES FOR USING THE MICROSCOPE

- 1) The lens is the most important component of the microscope. For this reason we advise using extreme care when handling the lens.
- 2) If the lens is dirty or dusty, the surfaces should be cleaned with an unsoiled cleaning paper or a soft cotton cloth. Never clean the glass surfaces of the lens directly with your fingers, dirty cloths, etc.
- 3) After use, the microscope should be kept covered at all times or returned to its packaging to protect it against dust.
- 4) The microscope should be stored in a dry location. Condensation on the light source can reduce the intensity of the light.
- 5) When the microscope is not in use for an extended period of time, the batteries for the light should be removed from the instrument.

INTRODUCTION TO THE WORLD OF THE MICROCOSMOS


There are a great number of living creatures in our world. Some are large and can be easily seen with the naked eye; others are so tiny that a million of them could easily fit on the head of a pin. We can observe such living creatures only with a microscope.

The microscope was invented many years ago, and since then research has opened an extraordinary universe that exceeds the bounds of our imagination with its fascinating beauty. Today, the microscope is used in nearly every scientific department, from the basic levels of biology to the highly specialised disciplines of astrophysics. The microscope provides students with a better understanding of the functions of complex life forms and organisms, as well as of the composition and structure of the materials that surround us. Your microscope will make your hobby more enjoyable, and it will provide you with an introduction to a more thorough study of a wide variety of sciences.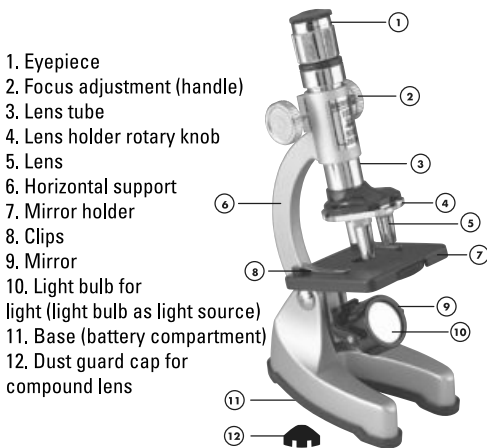


Operating Instruction

1. Tip the horizontal support and set the reflector so that the light is completely captured by the mirror.
2. Look through the eyepiece and check whether the light is completely reflected by the mirror; the microscope is now ready for use.
3. Place the previously prepared thin section onto the support plate and secure it with the retaining clips.
4. Now select a magnification. The magnification factor increases as the lens gets longer. Observation generally begins with the smallest magnification factor.
5. Turn the lens holder rotary knob for changing the magnification factor until you hear it click into position.
6. Turn the focus adjustment to move the lens down, getting it as close as possible to the section without touching it. At the same time, look through the eyepiece and adjust the focus adjustment in the opposite direction until the image comes into focus.
7. If the room in which the observation is taking place is weakly lit with sunlight or artificial light, or if the image cannot be clearly seen when you use a large magnification factor, the reflector should be turned over. The light bulb then lights up automatically to enable better observation.
8. The light source is a light bulb that is supplied with power from two 'AA' electric torch batteries (dry cells) on the back of the base.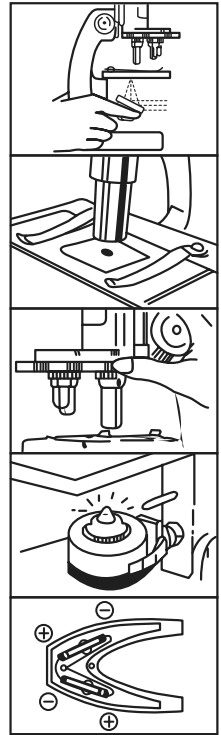

USE AS A PROJECTION DEVICE

Remove the eyepiece and tip the lens tube so that it is pointed at a white surface that is not more than a metre away. The best results are achieved in a darkened room.

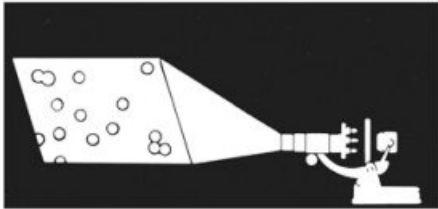ZOOM MODELS:

On the side opposite the lens tube, the numbers 12X and 18X appear.

Turn the silver-edged ring anti-clockwise to increase the magnification factor up to 18X.

Assuming you use the lens holder with magnification factor 12X and an eyepiece with magnification factor 12X, the optical system then has a magnification of $12 \times 12 = 144$. That means that the observed object is magnified 144 times.

If the zoom lens is turned around and set to the magnification factor of 18X, you obtain a combined magnification of $12 \times 18 = 216$.

ZOOM-LENS
12X-18X

CREATING SPECIMENS FOR OBSERVATION

If the sample piece is not thin and transparent, it cannot be observed under the microscope, since the light of the reflector and of the light source cannot penetrate it. Pollen fibres, wool or salt can be readily observed under the microscope and do not need to be covered with glass.

1) Provisional specimen

Clean the glass slide, cut off a thin slice from the sample with a razor blade or a similar tool (note: the edge is extremely sharp and needs to be handled with great caution) and place it onto the centre of the glass slide using tweezers. Place a bit of water onto the sample with a dissection needle (note: the dissection needle has a very sharp point and needs to be handled with great caution) and carefully cover the sample so that no air bubbles are enclosed with it. Remove any superfluous water using blotting paper. The sample can now be observed under the microscope.

USE OF THE THIN-LAYER SLICER

Place a sample that is to be cut into thin slices for observation into the opening of the thin-layer slicer (note: the instrument contains a very sharp edge and needs to be handled with great caution). Turn the blade. Thin slices of the sample material are created.

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(GB) Note on environmental protection:

After the implementation of the European Directive 2002/96/EU and 2006/66/EU in the national legal system, the following applies:

Electric and electronic devices as well as batteries must not be disposed of with household waste. Consumers are obliged by law to return electrical and electronic devices as well as batteries at the end of their service lives to the public collecting points set up for this purpose or point of sale. Details to this are defined by the national law of the respective country.

This symbol on the product, the instruction manual or the package indicates that a product is subject to these regulations.

By recycling, reusing the materials or other forms of utilising old devices/Batteries, you are making an important contribution to protecting our environment.

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Hama GmbH & Co KG